

## Technical Bulletin – Change to oxygen analyser replacement kit for NITROFLOW LAB Mobile Nitrogen Gas Generator

Dear colleagues, please be advised that the oxygen analyser replacement kit, part number 159.005574, will no longer include the housing assembly. The standard replacement kit will now contain the oxygen sensor unit only.

When replacing the oxygen sensor please follow the instructions below:

- 1) Take the analyser assembly and line up the screws by rotating the grey disk – Image A
- 2) Unscrew the 3 screws to free the sensor – Image B
- 3) Remove the grey housing – Image C
- 4) Gently pull the sensor away from the circuit board – Image D
- 5) Take the replacement sensor and use the guiding pins to plug the sensor to the circuit board – Image E
- 6) Once the new sensor is in place, reassemble the unit by aligning the 3 screws and tightening – Image F




*Image A*




*Image B*




*Image C*

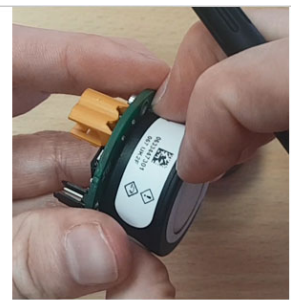

*Image D*

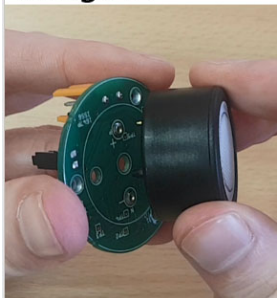

*Image E*

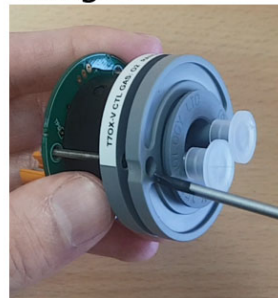

*Image F*

